


# Single-Walled Carbon Nanotube Dispersant

## Photochemically Controllable Dispersion and Aggregation

### Advantages

- Strong interaction to the surface of single-walled carbon nanotube (SWCNT)
- Positive charge on the stilbene core enables SWCNT to disperse in aqueous solution
- UV irradiation makes closure of the stilbene moiety and elimination from the SWCNT surface which enables SWCNT to aggregate
- Precisely controllable dispersion by photo irradiation
- After the aggregation, the dispersant can be easily removed by filtration, etc.

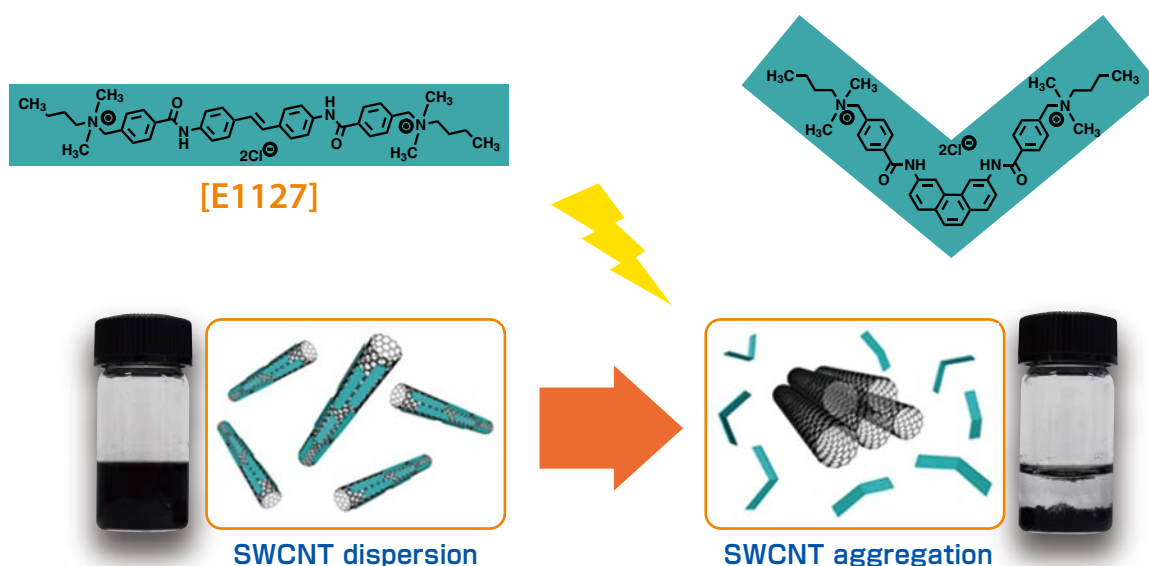

### Possible Application

- Purification of SWCNT by combining dispersion in solution, centrifugation and photo-induced aggregation
- Development of new composite material for CNT-based film and paint

References Y. Matsuzawa, M. Yoshida, H. Ohyama, H. Kato, Patent JP5552641.  
 Y. Matsuzawa, H. Kato, H. Ohyama, D. Nishide, H. Kataura, M. Yoshida, *Adv. Mater.* **2011**, 23, 3922.  
 Y. Matsuzawa, Y. Takada, T. Kodaira, H. Kihara, H. Kataura, M. Yoshida, *J. Phys. Chem. C* **2014**, 118, 5013.

**4,4'-[1,2-Ethenediylbis(4,1-phenyleneiminocarbonyl)]bis(N-butyl-N,N-dimethylbenzenemethanaminium) Dichloride**

1g [E1127]

# Single-Walled Carbon Nanotube Dispersant

## (1) Preparation of dispersions (Low power sonication)

3.6 mg of **E1127** was dissolved in 3 mL of pure water (18.8 M $\Omega$ -cm, TOC < 3 ppb) or D<sub>2</sub>O using a glass vial (ca. 0.12 wt%). The solution was then sonicated (80 W, 35 kHz) at room temperature. Heating by a dryer is also a useful method for solvation. 1.96 mg of SWCNTs was added into the solution of **E1127**. The **E1127**/SWCNT mixture was sonicated (80 W, 35 kHz, 1 h) at room temperature. The black-colored dispersion was then centrifuged to remove impurities such as amorphous carbons and metal particles (28500 x g, at 22 °C, 3 hrs). Upper 70 % of supernatant was corrected as a well-dispersed SWCNT dispersion.

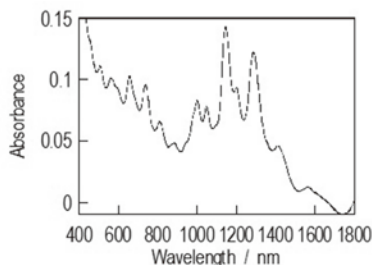


Figure 1. UV-vis-NIR absorption spectra of the **E1127**/SWCNT dispersion treated by low power sonication (in D<sub>2</sub>O, light pass length 0.5 mm)

## (2) Preparation of dispersions (High power sonication)

10 mg of **E1127** was dissolved in 20 mL of pure water (18.8 M $\Omega$ -cm, TOC < 3 ppb) or D<sub>2</sub>O using a glass vial (ca. 0.05 wt%). The solution was then sonicated (80 W, 35 kHz) at room temperature. Heating by a dryer is also a useful method for solvation. 7.0 mg of SWCNTs was added into the solution of **E1127**. The mixture was sonicated (80 W, 35 kHz, 1 h) at room temperature. Subsequently, the dispersion was subjected to ultrasonication (60 W, 19 kHz, 4 hrs) at room temperature. The black-colored dispersion was then centrifuged to remove impurities such as amorphous carbons and metal particles (28500 x g, at 22 °C, 3 hrs). Upper 70 % of supernatant was corrected as a well-dispersed SWCNT dispersion.

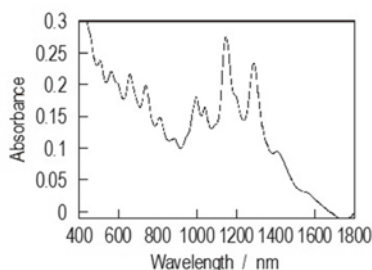


Figure 2. UV-vis-NIR absorption spectra of the **E1127**/SWCNT dispersion treated by high power sonication (in D<sub>2</sub>O, light pass length 0.5 mm)

## (3) Photoirradiation and precipitation

A well-dispersed SWCNT dispersion was transferred to a dialysis tube (cutoff MW of 1000) to remove excess amount of **E1127** from the dispersion. The tube filled with the dispersion of **E1127**/SWCNT system was put into a glass beaker with 3 L of pure water (18.8 M $\Omega$ -cm, TOC < 3 ppb) and was stirred overnight. The dialyzed dispersion was transferred into quartz cells and irradiated by an LED light source (365 nm, 100 mW/cm<sup>2</sup>) for several hours under magnetically stirring. After 3 hrs of light irradiation, precipitation of SWCNTs was completed for the dispersion treated with low power sonication. The SWCNT dispersion prepared using high power sonication took 7 hrs to recognize precipitation.

### Ordering and Customer Service

#### TCI AMERICA

Tel : 800-423-8616 / 503-283-1681  
Fax : 888-520-1075 / 503-283-1987  
E-mail : Sales-US@TCIchemicals.com

#### TCI EUROPE N.V.

Tel : +32 (0)3 735 07 00  
Fax : +32 (0)3 735 07 01  
E-mail : Sales-EU@TCIchemicals.com

#### TCI Deutschland GmbH

Tel : +49 (0)6196 64053-00  
Fax : +49 (0)6196 64053-01  
E-mail : Sales-DE@TCIchemicals.com

#### Tokyo Chemical Industry UK Ltd.

Tel : +44 (0)1865 784560  
Fax : +44 (0)1865 784561  
E-mail : Sales-UK@TCIchemicals.com

#### TCI Chemicals (India) Pvt. Ltd.

Tel : 1800 425 7889 / 044-2262 0909  
Fax : 044-2262 8902  
E-mail : Sales-IN@TCIchemicals.com

#### 梯希爱(上海)化成工业发展有限公司

Tel : 800-988-0390 / 021-67121386  
Fax : 021-6712-1385  
E-mail : Sales-CN@TCIchemicals.com

#### TOKYO CHEMICAL INDUSTRY CO., LTD.

Tel : +81 (0)3-5640-8878  
Fax : +81 (0)3-5640-8902  
E-mail : globalbusiness@TCIchemicals.com

Availability, price or specification of the listed products are subject to change without prior notice. Reproduction forbidden without the prior written consent of Tokyo Chemical Industry Co., Ltd.